

Mounting a Coolpix 5000 onto a Microscope Reference Sheet*

Please refer to the Coolpix 5000 User's Manual for page numbers indicated below.

Mounting the Coolpix 5000 to a microscope trinocular photo tube or beamsplitter

1. With the camera power off, open the card-slot cover and make sure a CompactFlash™ Card or IBM Microdrive® is properly inserted into the camera. **See p.36**
2. Slowly thread the UR-E6 Step Down Adapter to the camera.
3. Carefully thread the MDC-A Relay Lens onto the UR-E6 Step Down Adapter. Likewise, thread the C-mount onto the MDC-Relay Lens securely.

Note that over-tightening the adapter may damage threads of both adapter and apparatus.

4. Attach the C-mount adapter onto the microscope's trinocular or beamsplitter port. Secure the C-mount into place.
5. Adjust the LCD screen to the desired viewing position and angle.

Note: Use a 1x C-mount (e.g. Nikon ISO 1x C-mount, #83043) to ensure the maximum field of view; EH-52 AC Adapter is recommended to eliminate battery consumption.



Mounting the Coolpix 5000 into a 30mm eyepiece tube

1. With the camera power off, open the card-slot cover and make sure a CompactFlash™ Card or IBM Microdrive® is properly inserted into the camera.

See p.36

2. Slowly thread the UR-E6 Step Down Adapter to the camera.
3. Loosen the three setscrews of outer sleeve on the MDC-A Relay Lens and remove the Cover Ring.
4. Carefully thread the MDC-A relay lens to the UR-E6 Step Down Adapter.
Note that over-tightening the adapter may damage threads of both adapter and apparatus.
5. Remove an eyepiece from the microscope and insert the available end of the MDC-A Relay Lens adapter as replacement.
6. Adjust the LCD screen to the desired position and angle.
7. The weight of the camera may cause its positioning to become unstable.

UR-E6 Step Down Adapter



Cover Ring

Setscrew (x3)



Note: Nikon Eclipse and Nikon Stereomicroscopes are 30mm in diameter. Other Nikon models & microscopes from other manufacturers may vary. The EH-52 AC Adapter is recommended to eliminate battery consumption.

* This reference sheet is to be used in addition to the Coolpix 5000 User's Manual

Coolpix 5000 Camera Settings for Microscopy

1. Turn the camera on and select the infinity mode by pressing the *Manual Focus (MF)* button until the icon with two mountain peaks is displayed in the upper-right of the monitor display. Note that the flash turns off automatically when focus is set to infinity **See p. 77**
2. Hold down the *MODE* button and simultaneously turn the *COMMAND DIAL* to select **A**, *Aperture-Priority Auto* mode. The camera will set shutter speed automatically for best results.
3. Rotate the *COMMAND DIAL* to set the **F** value to the lowest possible f-stop number. The approximate range value (F2.5 - F4.0) is dependent on the zoom setting.
4. Use the zoom buttons (**W-T**) to fill the field of view. Use the LCD screen to determine best image. An image without black cut-off corners or circular images is generally possible at mid-range of the zoom setting.

Optional Settings in Menu Selection

Above the LCD display, press center-button above *Menu* for the Shooting Menu for camera settings available **only** in the **P**, **S**, **A**, and **M** mode and **only** when the camera is in shooting mode. The following are recommended options for photomicrography:

- ? **White Balance** The default setting is *Auto*. Select a white balance setting that matches the microscope light source; incandescent for tungsten or halogen illumination. Select *White Bal Preset* for unusual lighting conditions. **See p. 102-104**
- ? **Image Quality:** Image quality settings other than *Hi* are compressed as they are saved into memory. Compression selectively reduces the quality of an image. Certain illumination techniques, such as DIC, phase contrast, OCC, etc., may require special brightness and control adjustments for best results. **See p. 72-73**
- ? **Metering:** The default setting is *Matrix*. Setting it to *Spot* on a targeted sample with a very dark or a very bright background would ensure most appropriate exposure. **See p. 105**
- ? **Continuous:** The default setting is **S**, single shutter release. Using other settings, continuous sequence of pictures may be taken while the shutter-release button is held down. **See p. 108-109**
- ? **Saturation Control:** The *Black&White* setting will enhance black and white samples. **See p. 115**

